



Comprehensive mobile application to monitor your Coriolis Compact

- Remotely control your Coriolis Compact
- Get access to all functionalities in one mobile app
- Easily set up your collection protocol
- Available in English & French

CORIOLIS CONNECT

Coriolis Connect is a mobile app that can be used to control your Coriolis Compact instrument remotely. As soon as you start your Coriolis Compact air sampler, the instrument can be automatically detected by the app via Bluetooth. With the app, you can control your Coriolis Compact and easily register new protocols. During sampling, you can have access to a notepad section to input all the remarks of your protocol.

Whether working in harsh environments or in clean rooms, you will be able to step out and remotely control the Coriolis through the Coriolis Connect app without interfering with workers or the sampling environment.

Download all your sampling history, maximize your productivity, and complete your LIMS (Lab Information Management Systems) or ELN (Electronic Lab Notebooks).

Quick User Guide



1. Download the application on the website



2. Activate the bluetooth on your Coriolis and select it in the application




3. Edit the parameters



4. Start the cycle

Coriolis Connect is compatible with the Coriolis Compact firmware version v2.3.6 or latest version. Coriolis Connect is only available on Android devices.

Discover our comprehensive range of solutions

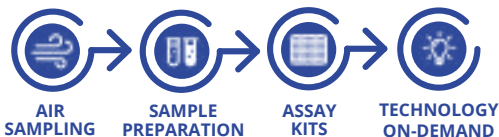

Bertin Technologies SAS
Parc d'activités du Pas du Lac
10 bis, avenue Ampère
78180 Montigny-le-Bretonneux, FRANCE
01 39 30 60 00

WWW.BERTIN-TECHNOLOGIES.COM